

Gly medium recipe

Stefano Varricchio - EPFL iGEM 2010

July 14, 2010

Where On the wetbench at the right of the fume hood

		Quantity	What
Material needed	For 1 liter:	25 g	Glycerol
		10 g	Yeast extract
		1 L	diH ₂ O
		1	Beaker

Steps

1. mix glycerol, Bacto yeast extract and dH₂O.
2. adjust pH until pH = 5
3. Sterilize by autoclaving.

Warnings

- Before doing a lot of bottles of medium try when they fit in the autoclave
 - Don't close the lid hermetically for autoclaving, the bottle will *explode*. Let air flow in and out the bottle.
 - Bacto yeast stinks a lot.
-